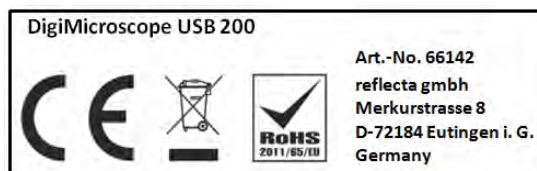
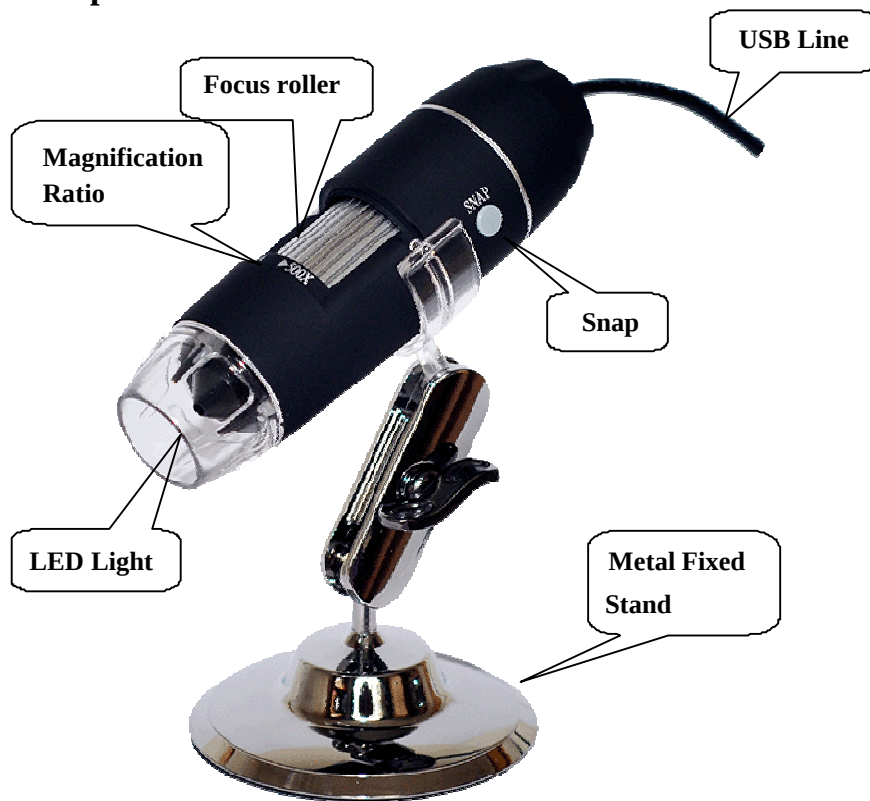
reflecta DigiMicroscope USB 200

User Manual

www.reflecta.de

Parts of Microscope

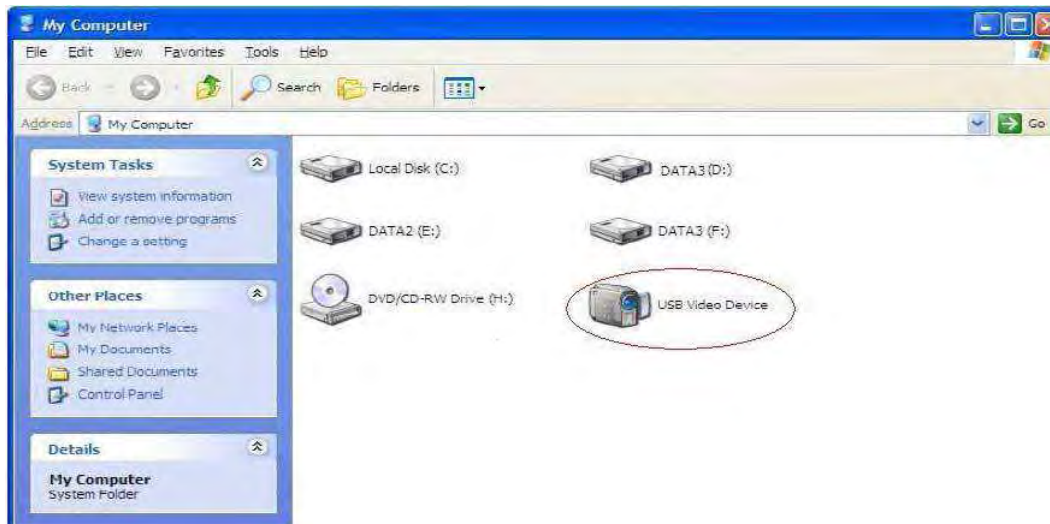
Notes before use

1. Don't disassemble the digital Microscope or change the interior parts, it can cause damage.
2. Don't clean the Microscope with alcohol organic solvents
3. Don't touch the lens with your fingers.
4. Avoid outdoor use if possible.
5. Storage temperature , 0° C ~ 40° C, Humidity:45%RH ~ 85RH%.
6. In case the product gets wet, leave PC connection immediately. And do not disassemble or dry by hair dryer. Send to repair center if the digit microscope was effected by liquid or other elements.
7. Measurement Data only for reference.

Scope of delivery

- Digital microscope with USB interface
- Stand
- Software CD
- Quick Start Guide
- Calibration ruler

Hardware system requirements

Windows XP, VISTA, WIN7, WIN8, WIN10 Pentium 1G, Celeron, AMD 1G & above,
128MB Memory, 150MB Hard Disc memory space, 16-bit & above
VGA, CD-ROM, USB2.0 or USB1.1.

How to install the Driver

The Digital Microscope is a free Driver Product, it can be worked immediately after connected to the computer with operation system above win2000 version.

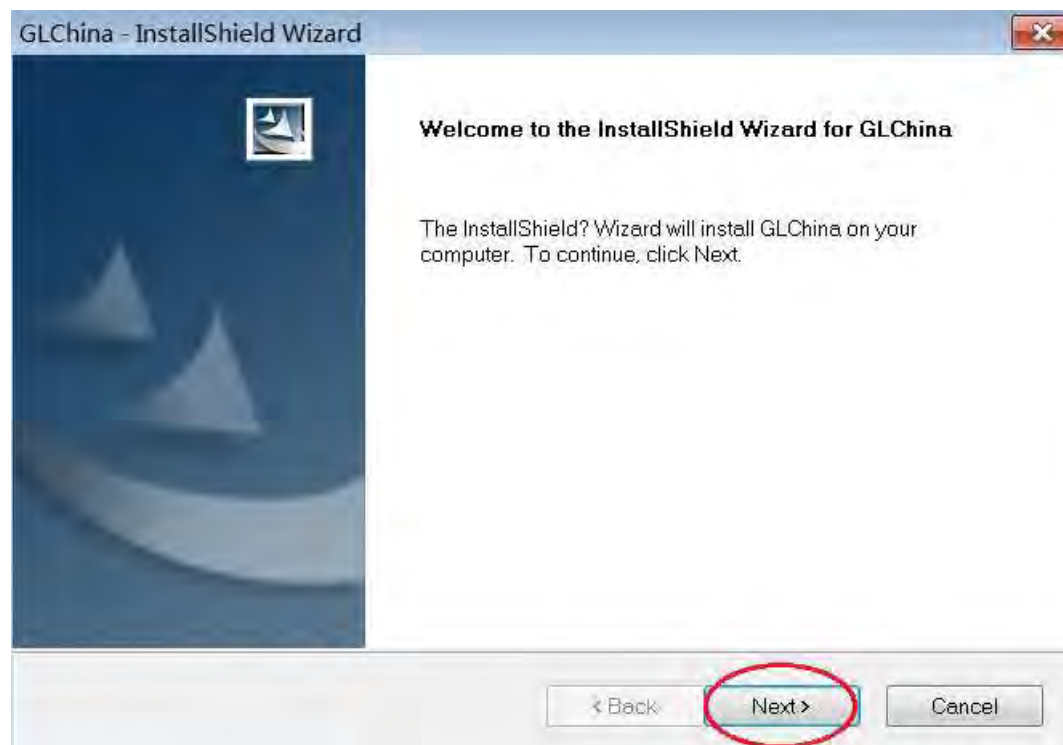
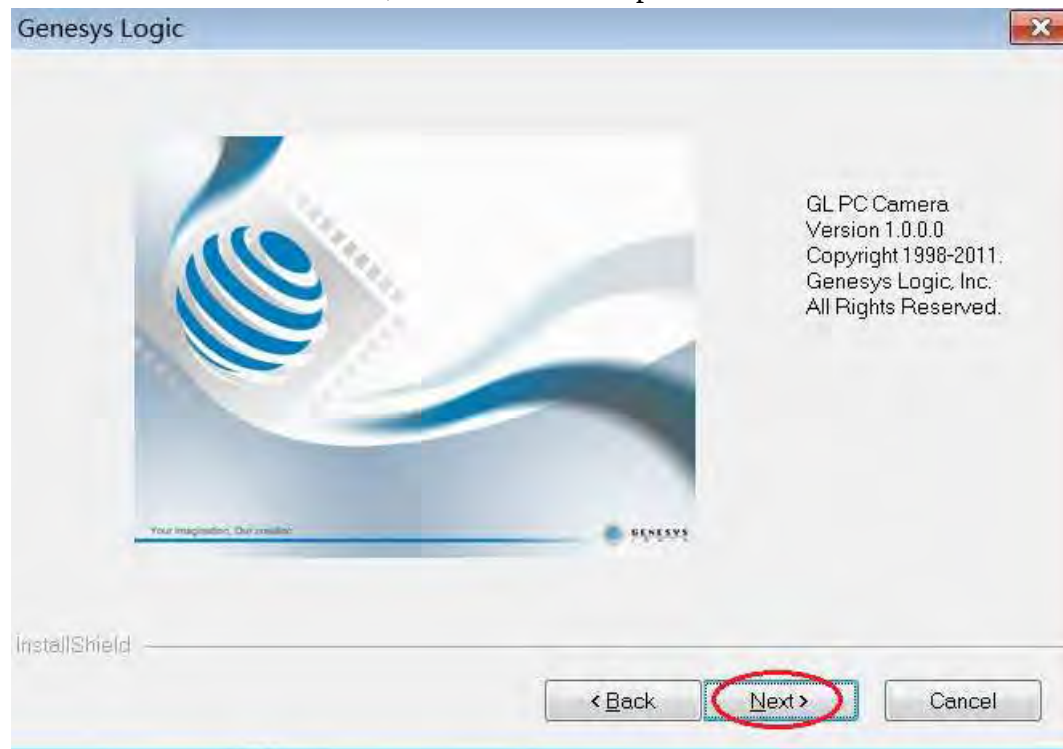
1. Connect the Portable digital microscope by USB2.0 cable to your computer, double click “my computer” icon on your computer desktop, below interface will be shown up.
2. Double click the Video device icon as above red marked in your computer. Now it is ready to use.

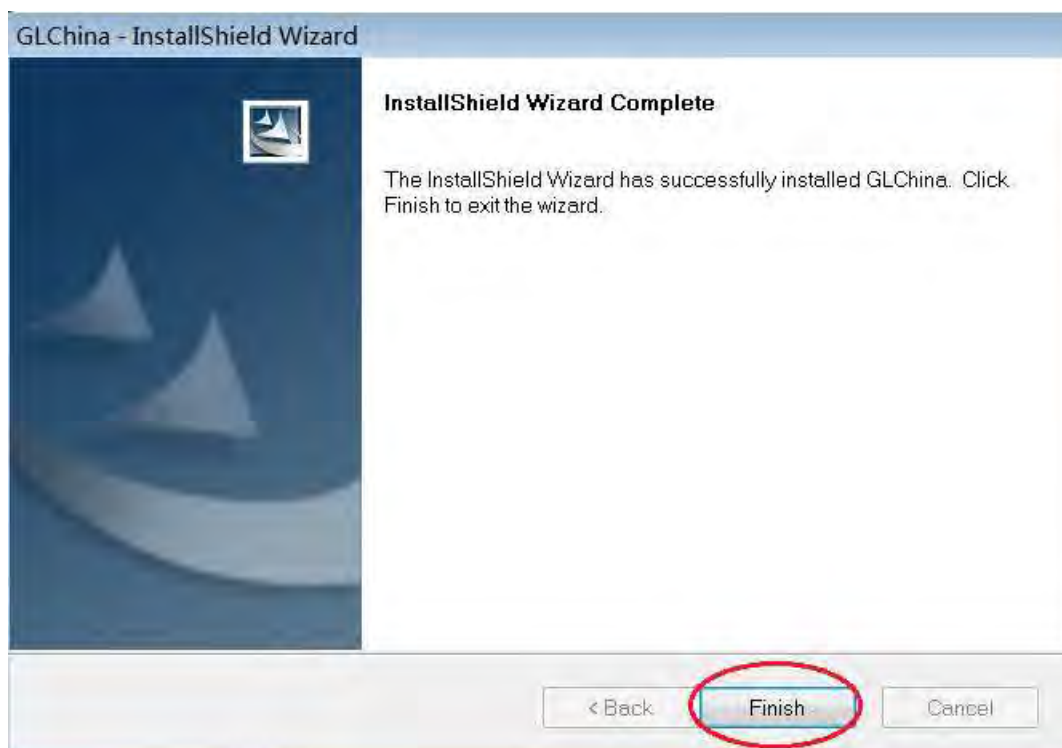
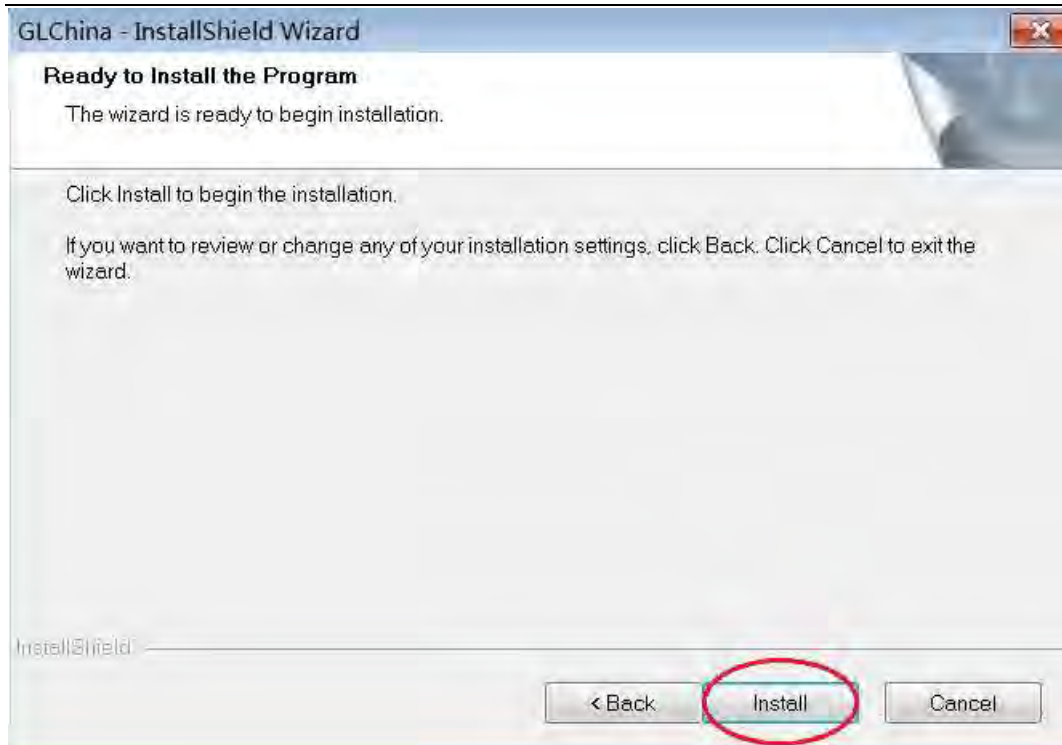
If you can't find the Video device, or if your computer can't recognize it,
Please follow below steps to install the driver.

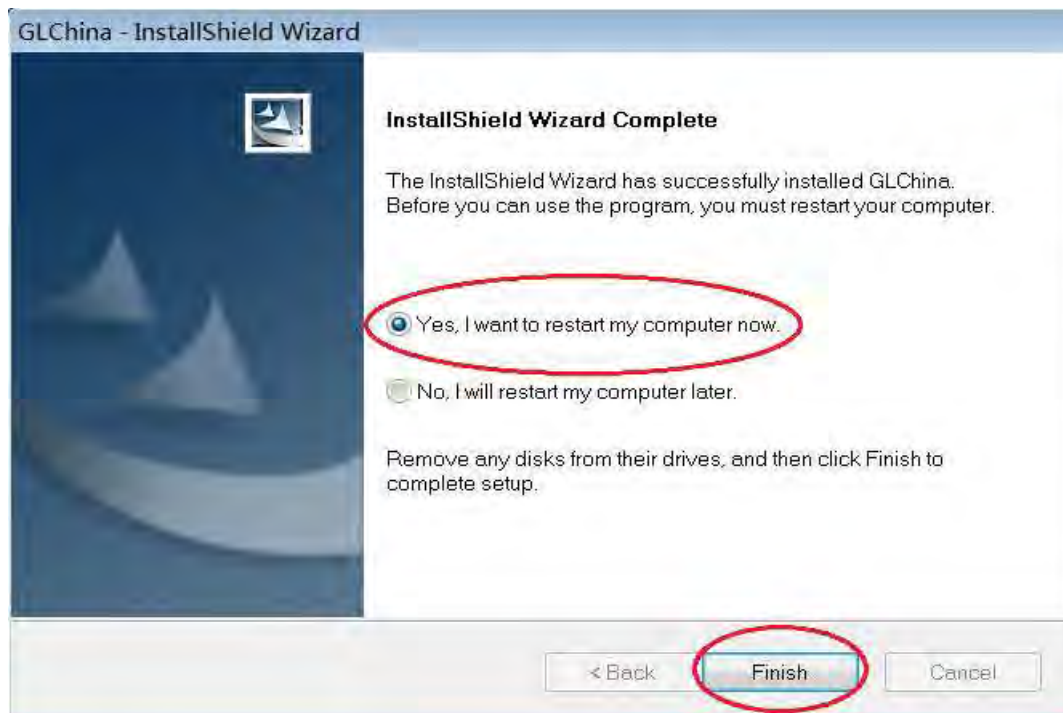
Insert the attached CD, open “My Computer”, and double click the “DVD/CD Driver” to open the CD.

Driver installation

Choose and install the Driver, then reset the computer

Reset the computer after driver installation at once;

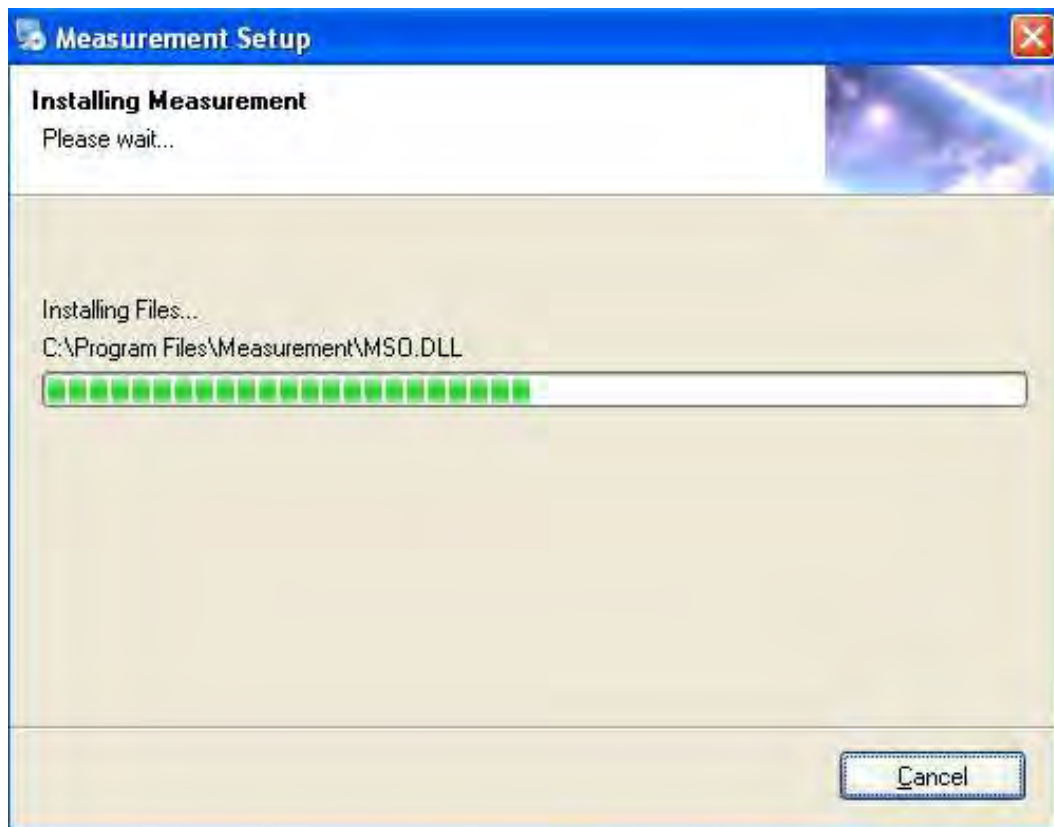
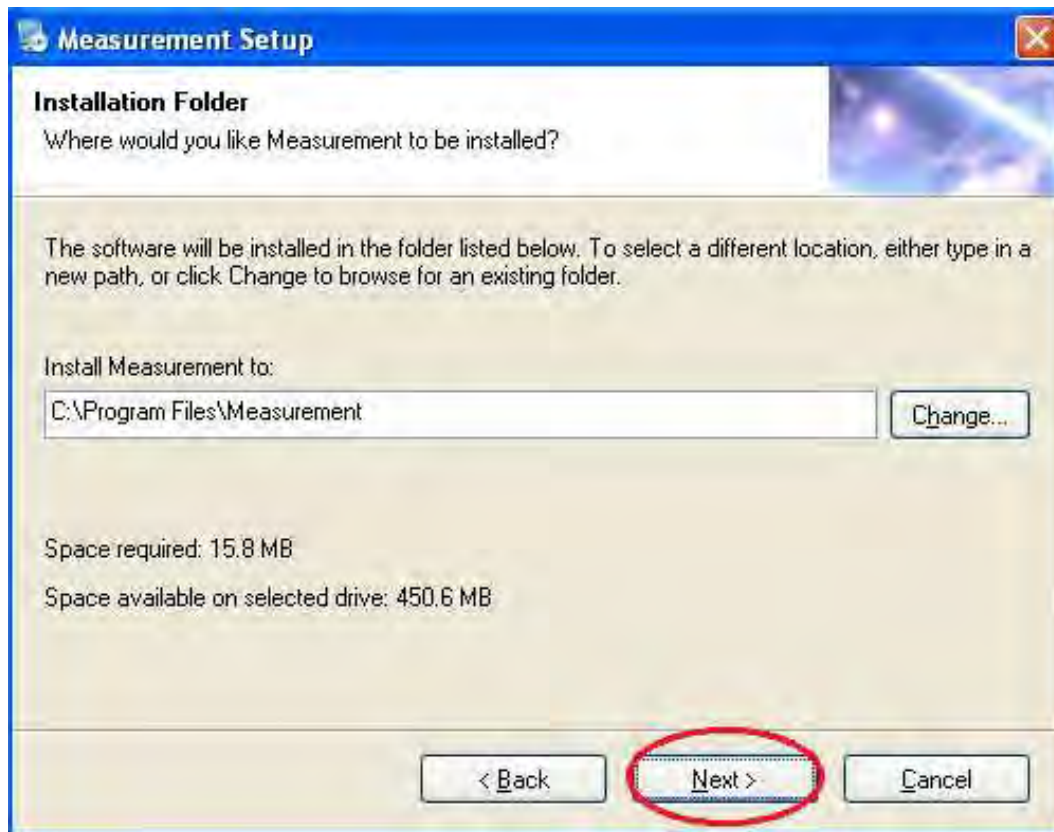
Installation may be blocked by the antivirus software, so let the antivirus software allow all processes, or close the antivirus before installing. If installation succeeds, you can see the icon  on the Toolbar, also the software desktop icon, double click to run it.

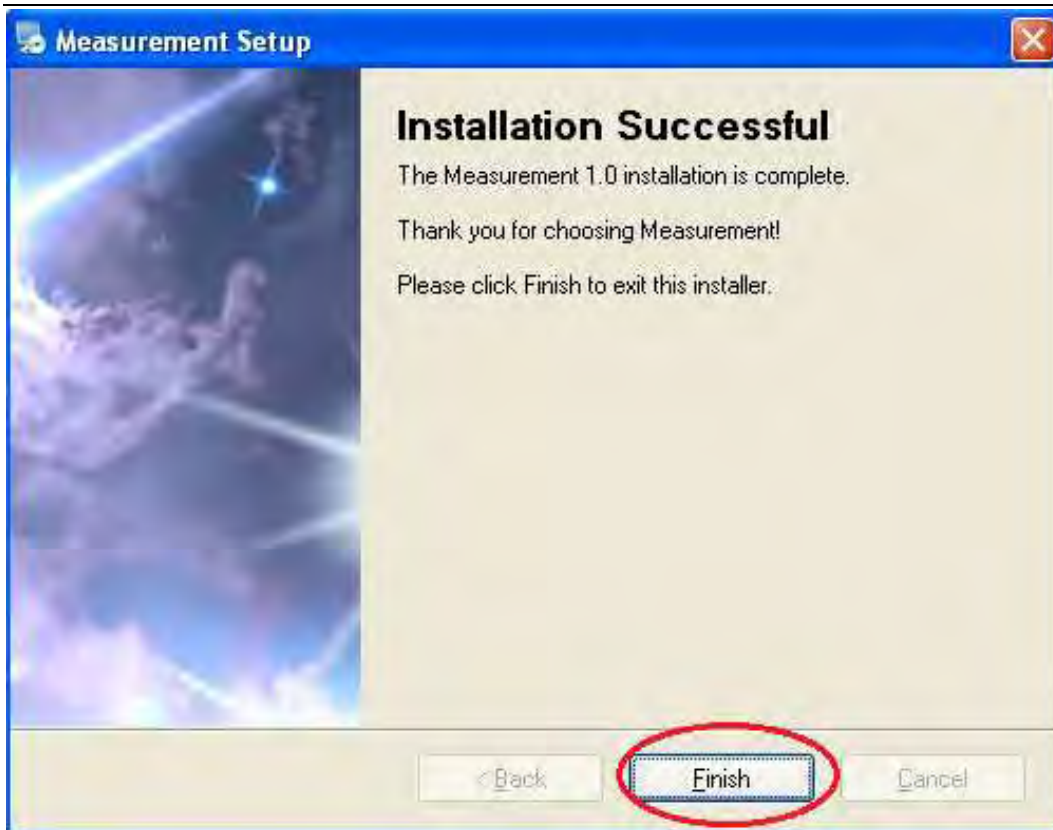
Now you can use the microscope. Long press the SNAP button on the microscope to take images, and the ZOOM button to have digital zoom function.

Measurement installation

Choose and install the Measurement



Click "Finish" and the installation completes.

Microsoft .NET Framework installation

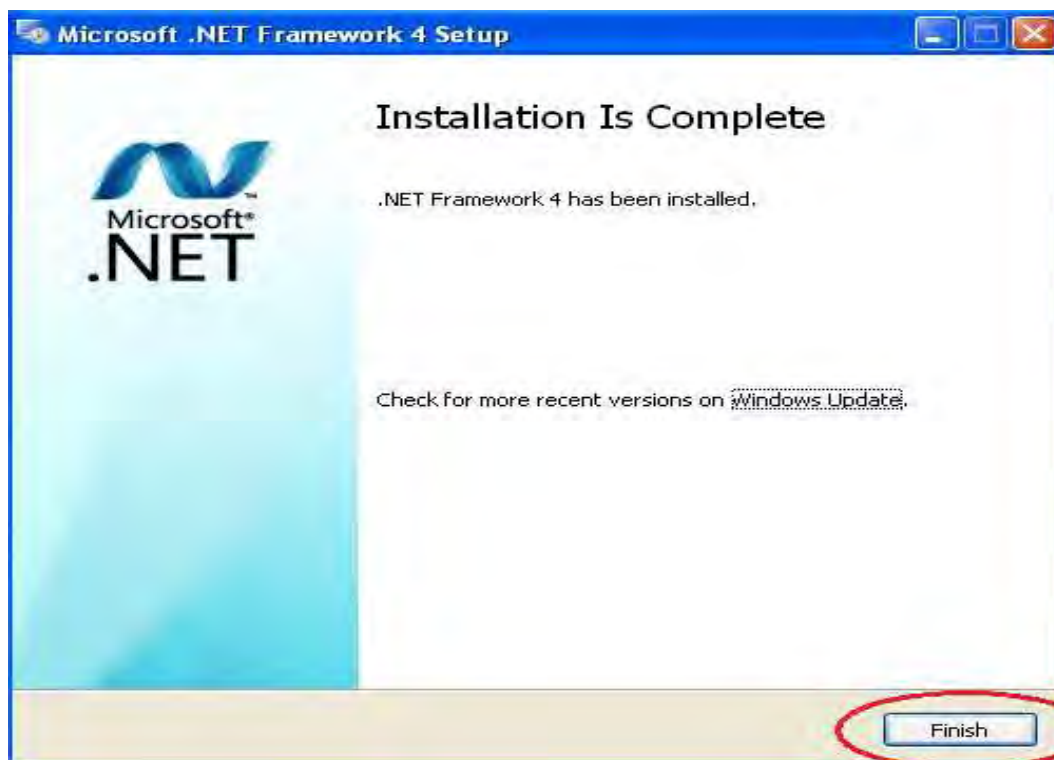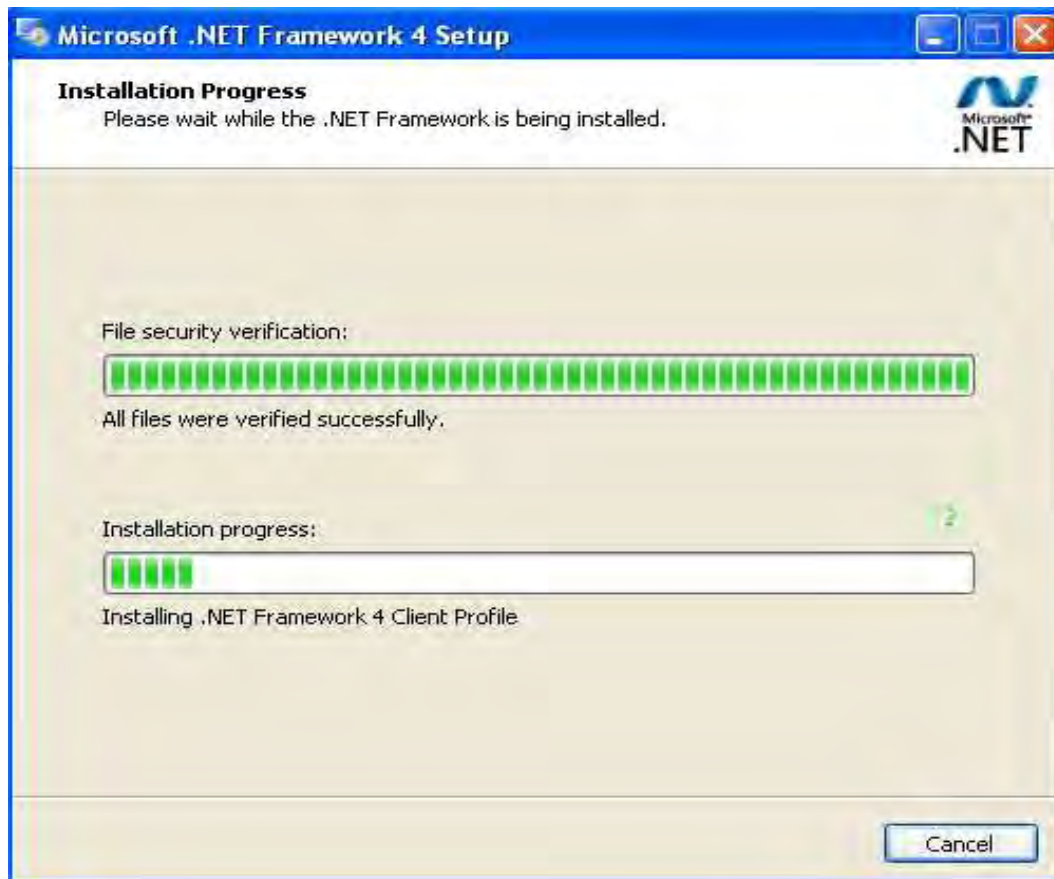If the PC is Windows XP or older version, .Net Framework 4.0 is needed. Install it.



After the installation completes, double click  at computer desktop to start using the software.

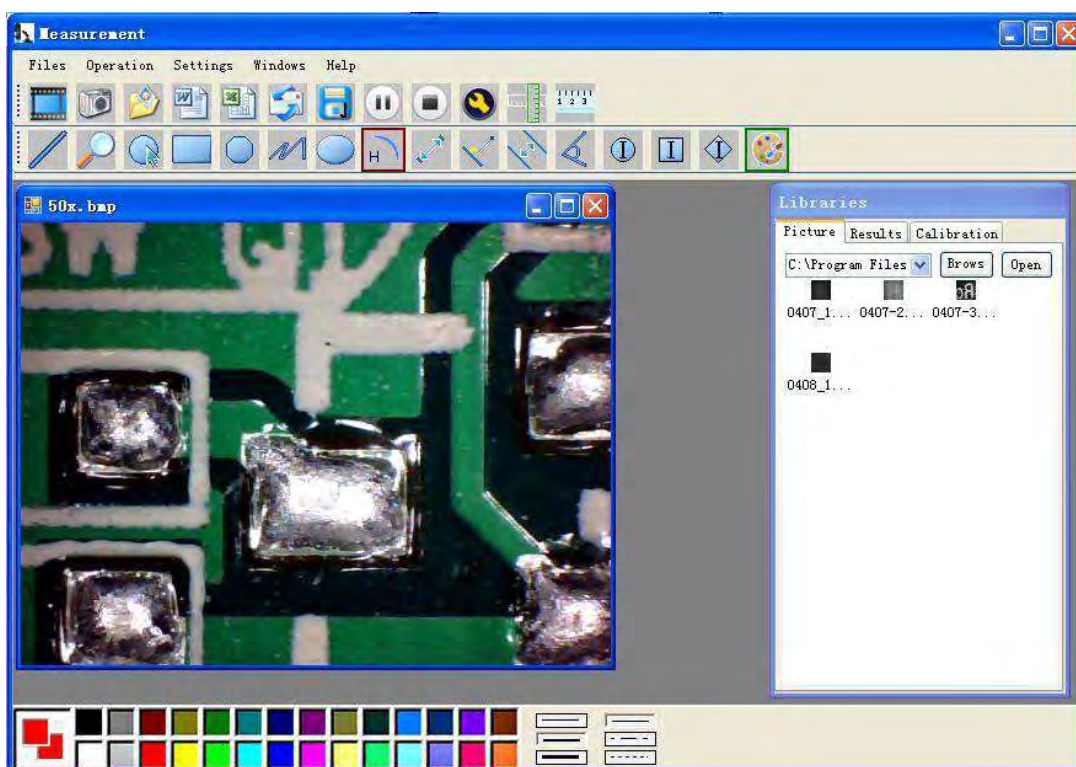Plug-in the devices, Display in operation window


Click  open the operation system

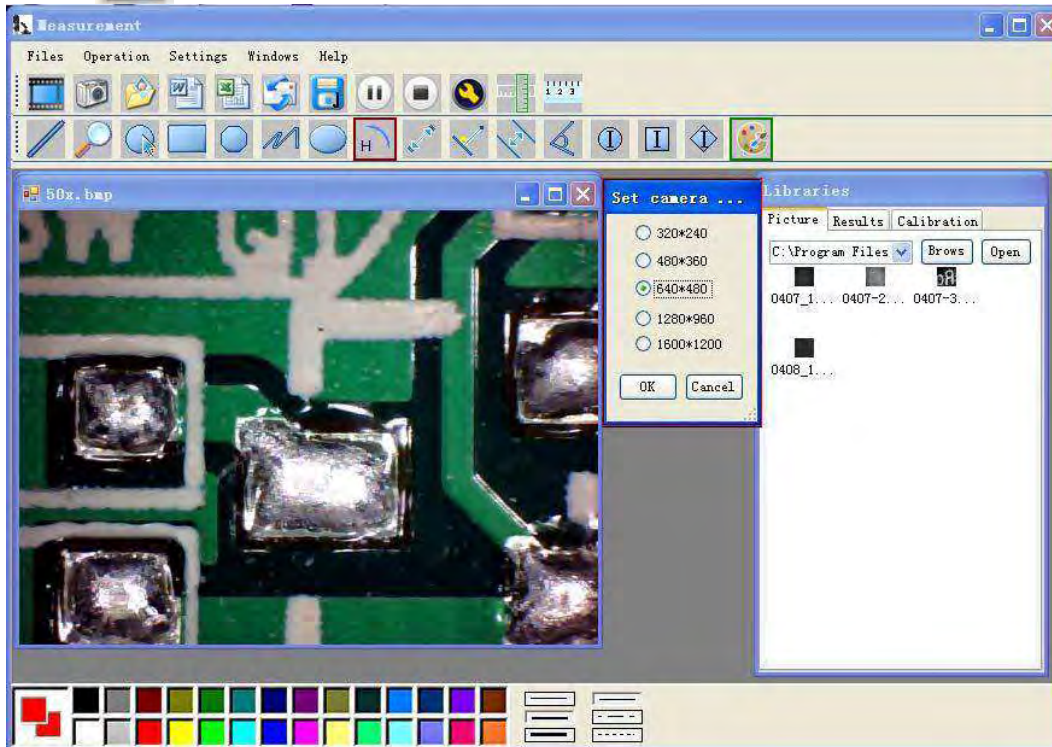
Choose the Image resolution from 320*240 to 1600*1200, default at VGA Mode 640*480

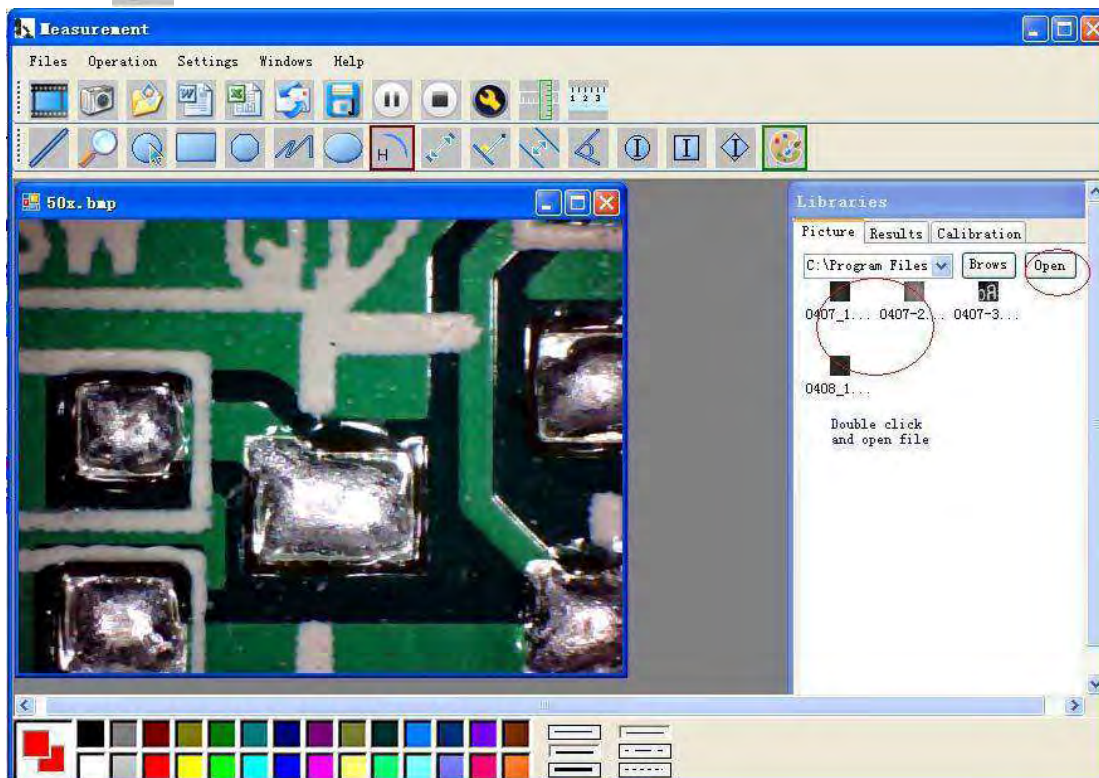Click  To use measurement function.

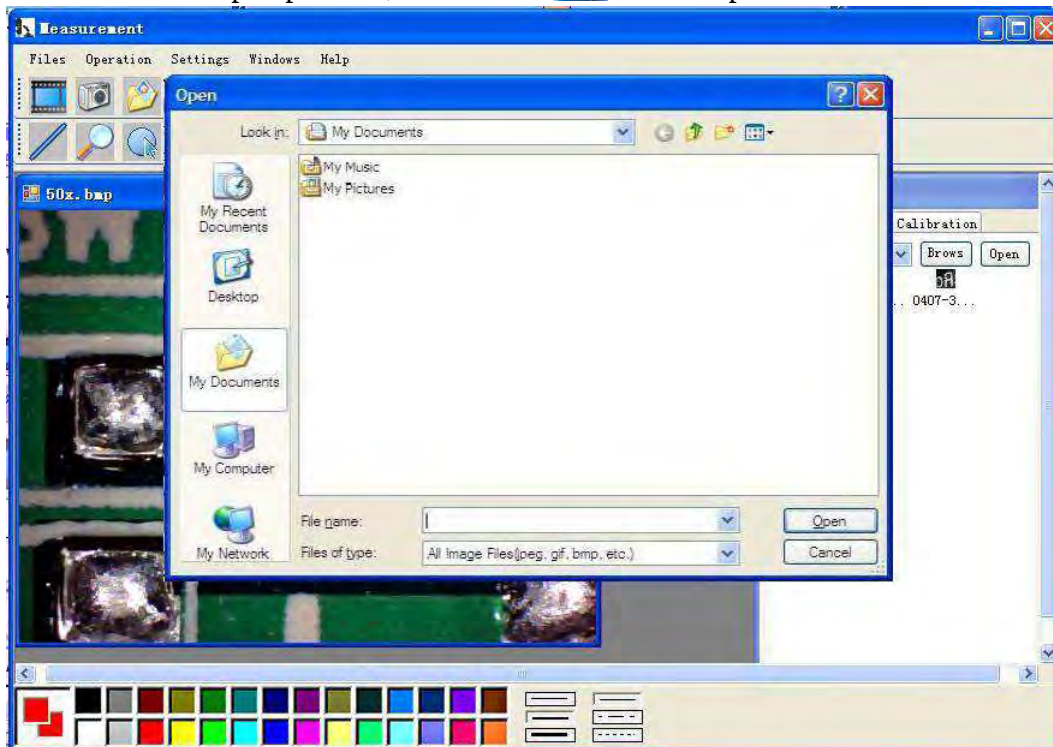
Click  to open pictures, and click  to save pictures

Video PLAY (Pause) ,STOP

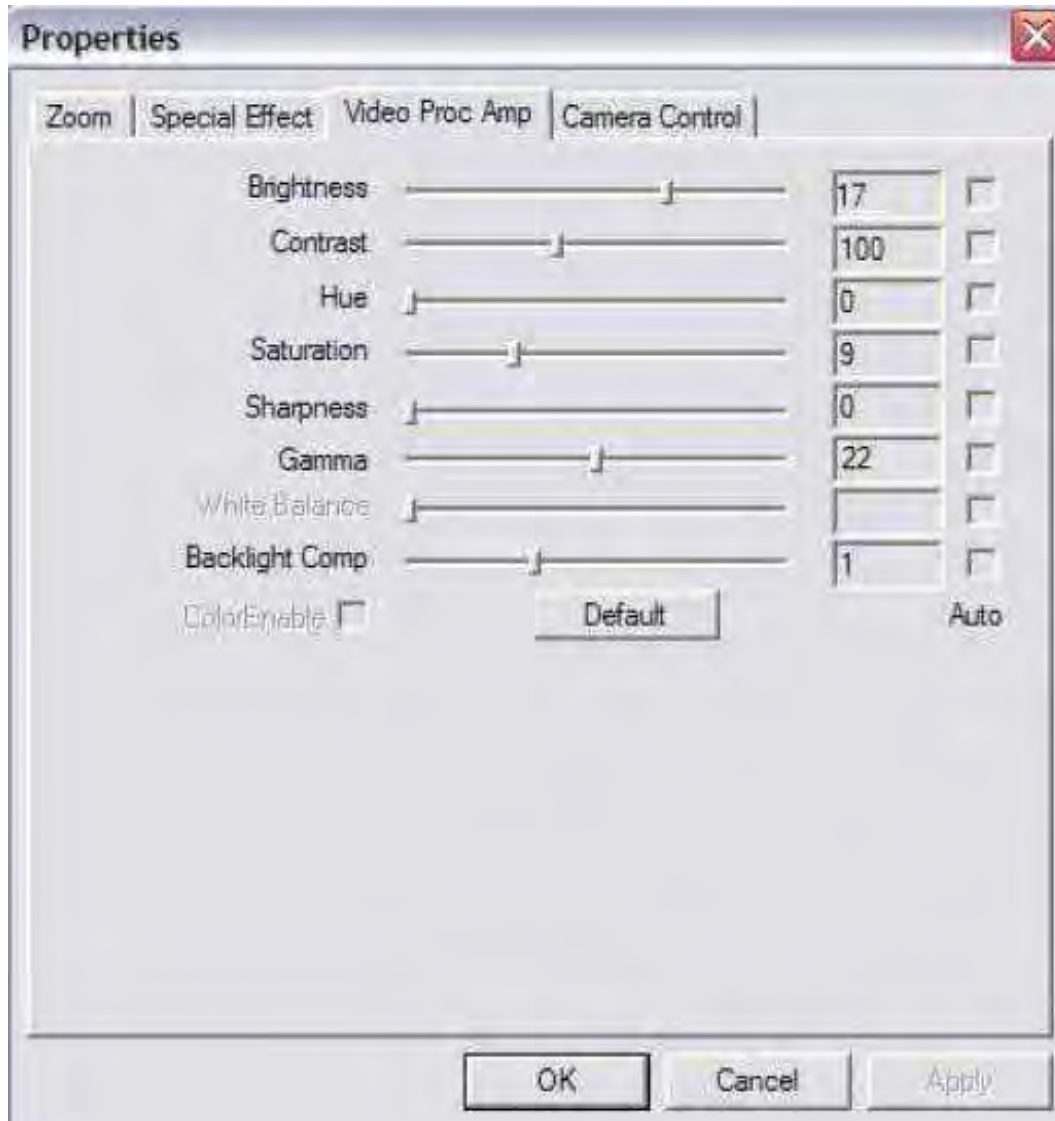
Image Parameter including Image setting and special effects

Magnifier

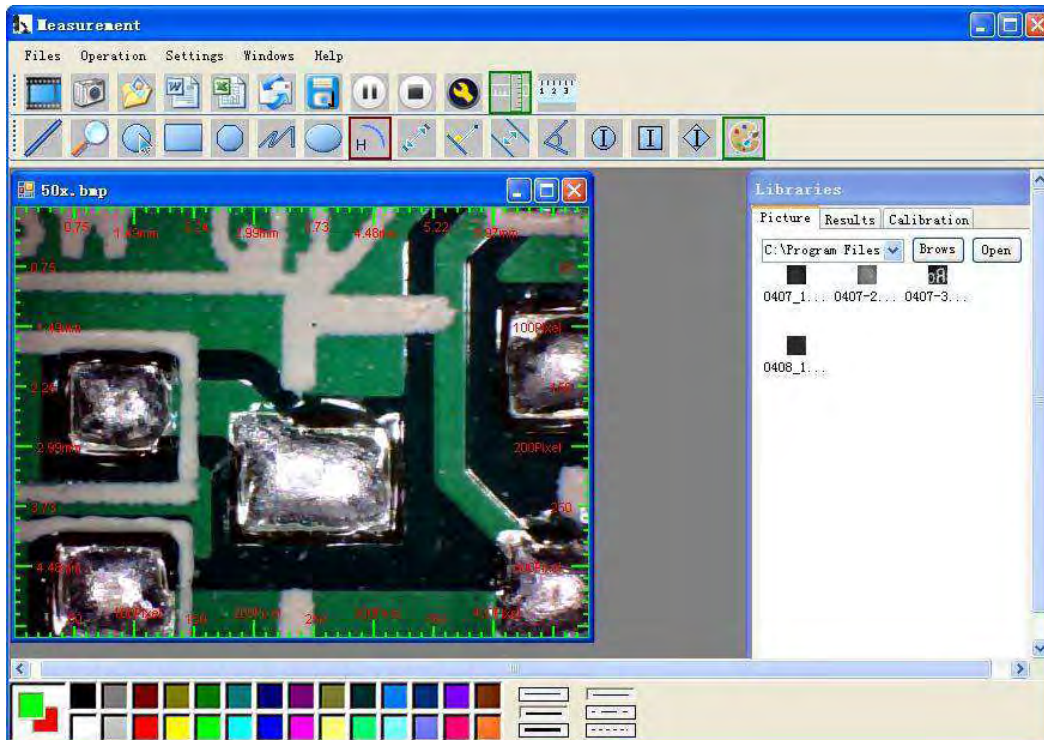
Click Ruler, as followings


Click start calibration make sure certain accuracy.

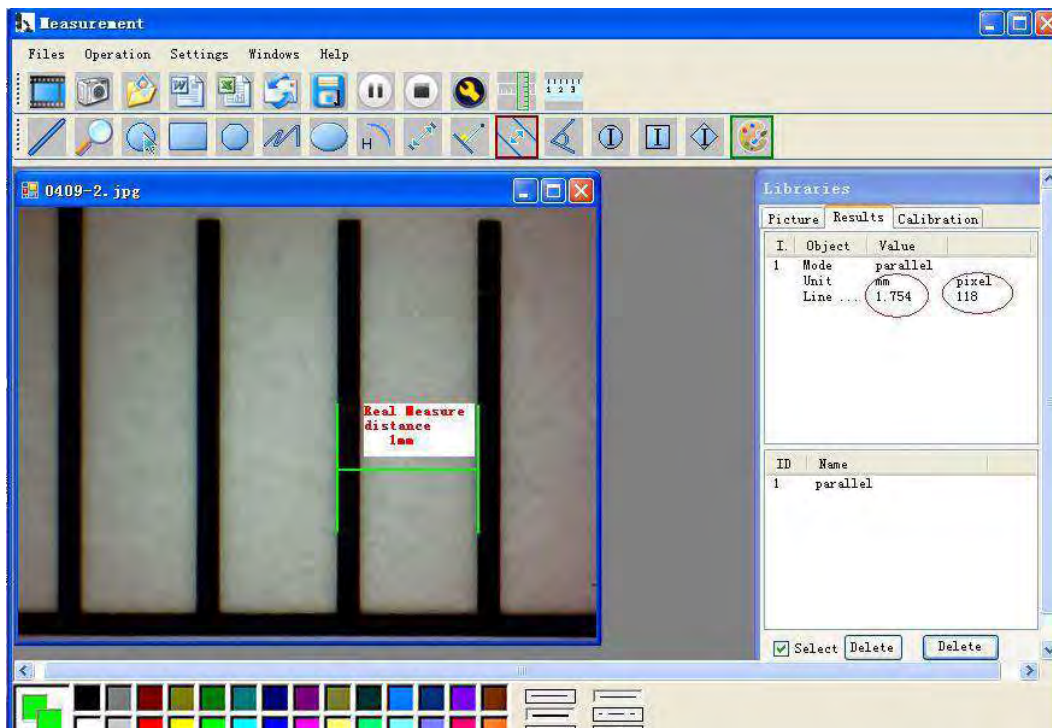Our experience told us that measurement could be controlled at 99% high precision. Measure tolerance could be caused from human eye distortion and defocus. More practice is helpful to measurements.

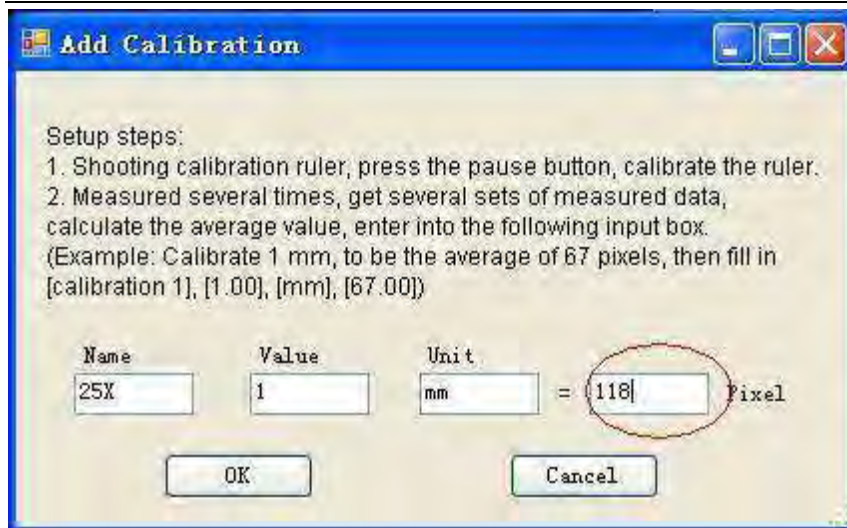
Key in the Measure data (Pixel), Unit, and Click' OK'

Warning: Please do calibration again once Magnification Ratio changed.

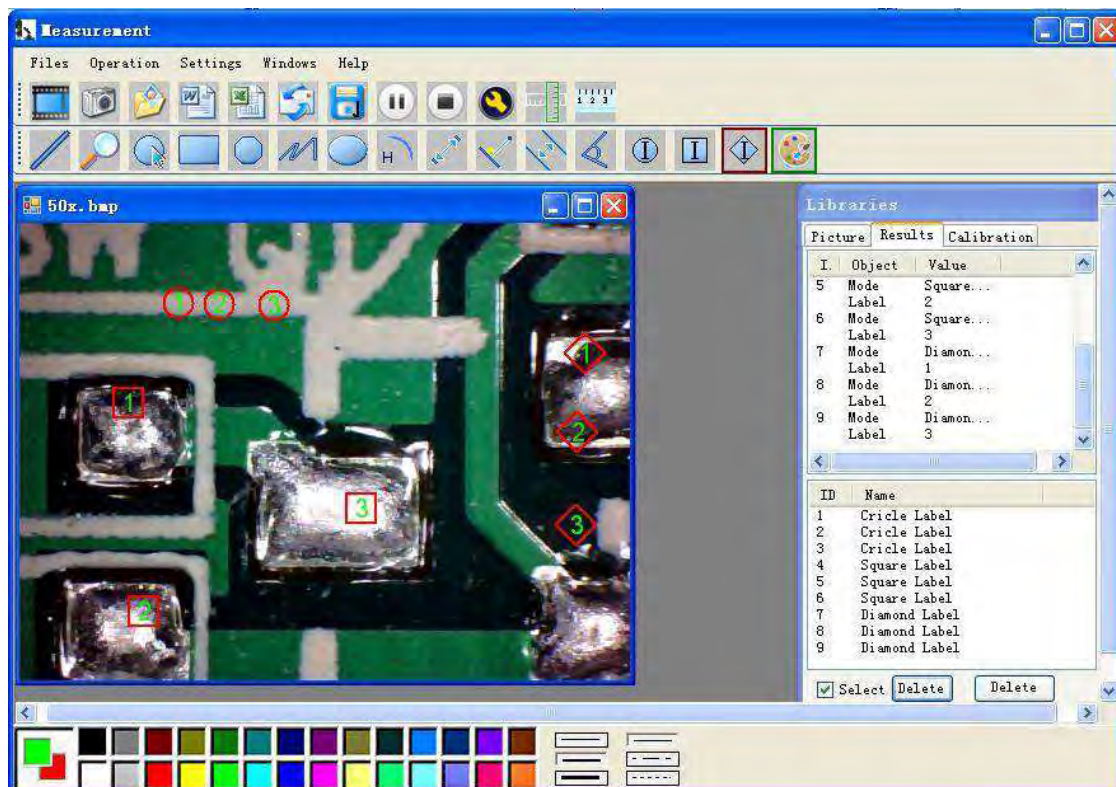Color setting of display character

Click to export word or excel format.

Example: Measure data under word file

Result output

Time: 2010-04-09 17:27

Data:

Mode	Unit	Line Length
parallel	mm	0.688
parallel	mm	非数字
parallel	mm	0.000

Mode	Unit	Angle
Angle	Degree	129.289
Angle	Degree	128.425
Angle	Degree	135.000

Picture:

From Left to right:

Line, ring, Square, Polygon, Set-line Length, Oval, Arc, Point to Point distance, Point to Line distance, Parallel line, Angle

Measurement

Files Operation Settings Windows Help

50x. bmp

Radius 0.696mm
Radiant 0.927mm

3.821 mm

1.115 mm

Libraries


I	Object	Value
8	Mode	Square...
	Label	4
9	Mode	Diamond...
	Label	4
10	Mode	Ellipse
	Unit	mm pixel
	Longe...	1.164 78
	Short...	0.567 38
	Perim...	2.720 182
	Area	0.519 2327

ID	Name
2	Cyclometry
3	Rectangle
4	Polyline
5	Arc
6	parallel
7	Circle Label
8	Square Label
9	Diamond Label
10	Ellipse

Select Delete Delete


For example 1

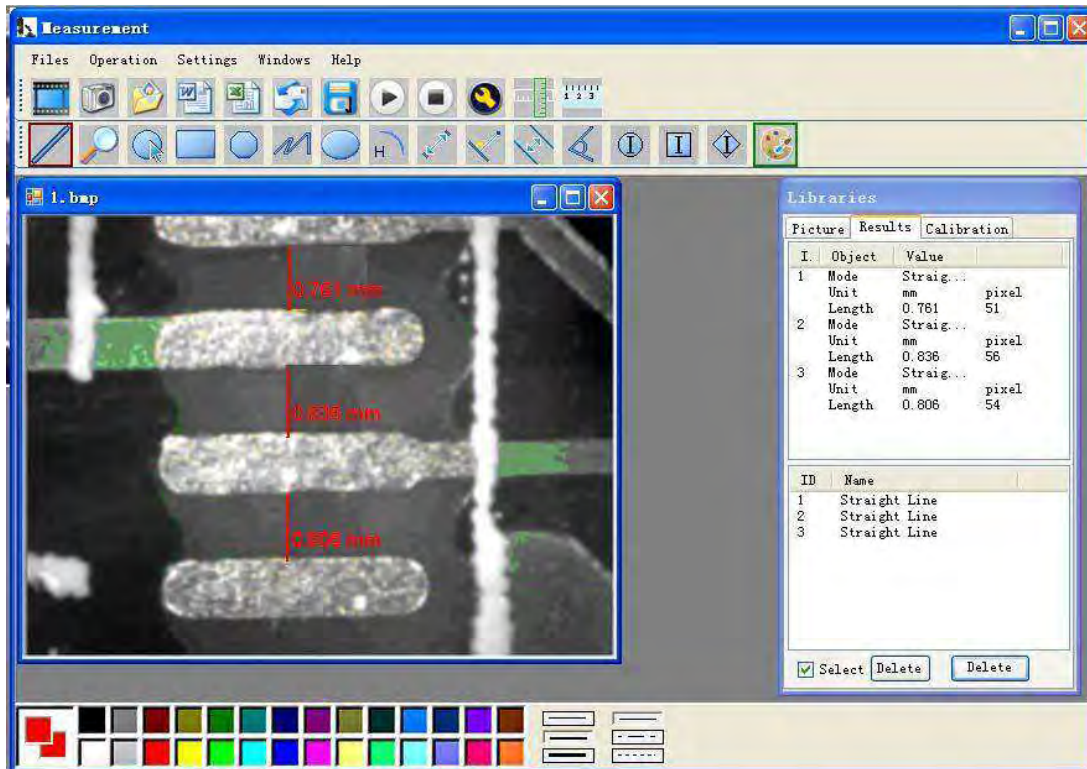Line measurement:


For example 2

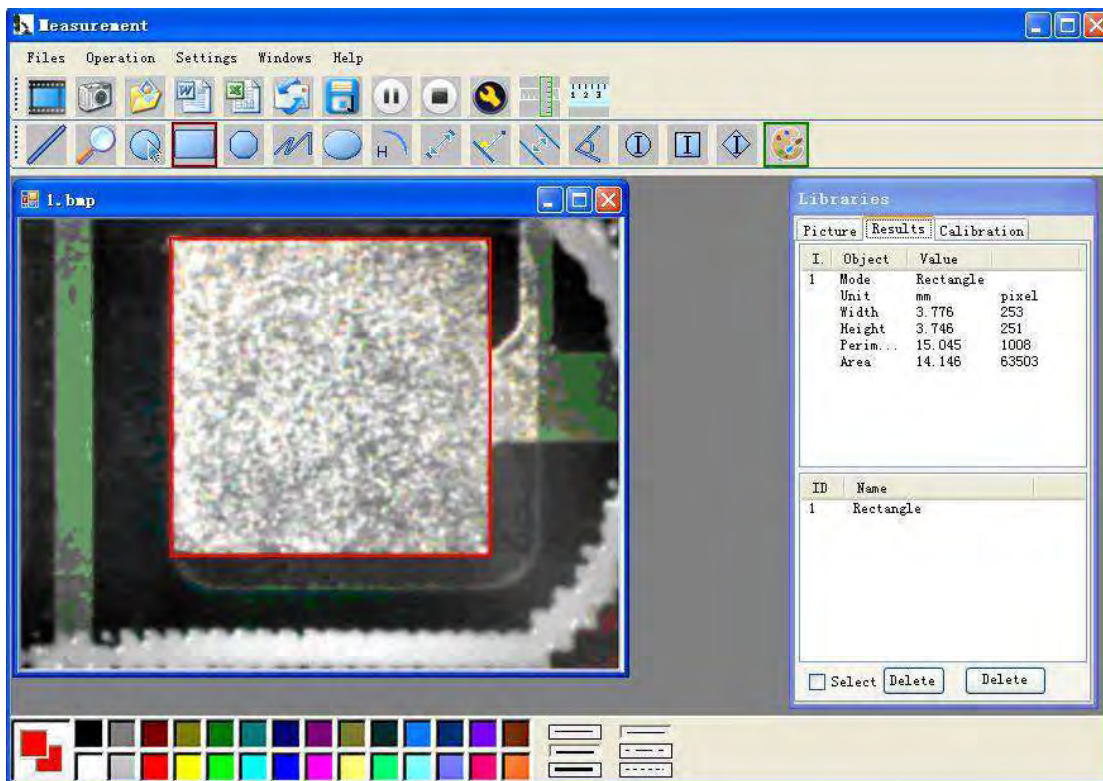
Rectangular measurements:

For example3
Ring measurements:

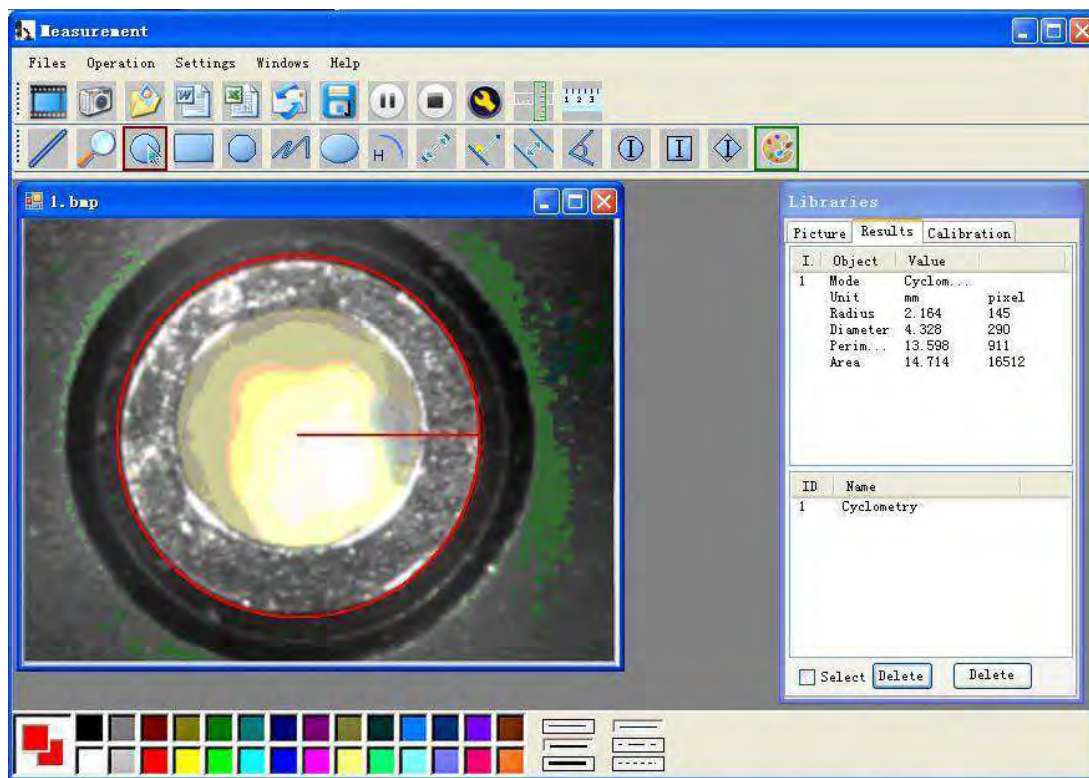For example 4
Arc measurements:

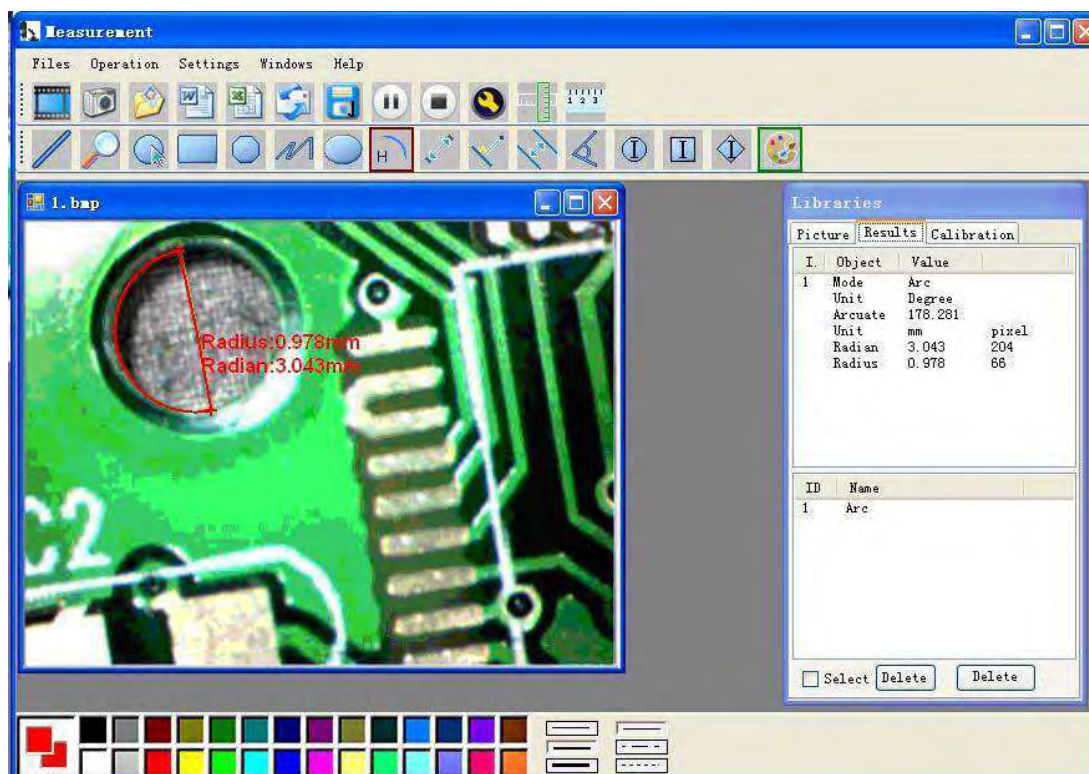This symbol on the product or in the instructions means that your electrical and electronic equipment should be disposed at the end of its life separately from your household waste.

There are separate collection systems for recycling in the EU.

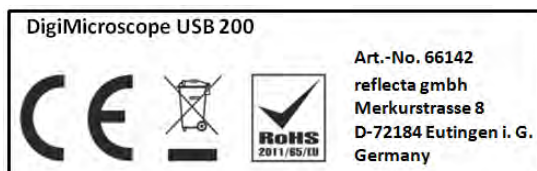
For more information, please contact the local authority or your retailer where you purchased the product.

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